

Elite™ Firefly Luciferase Assay Kit

CATALOG NUMBER: CA-L165, CA-L165-10

Updated on January 26, 2023

Description

Provided in the sizes of 10 mL (CA-L165, for one plate) and 100 mL (CA-L165-10, for 10 plates)

Application

- For firefly luciferase reporter assay

Features

- **Greater linear range**
- **Accurate**
- **Robust:** Excellent signal to noise (basal) ratio. Stable assay signal.
- **Easy to use:** Homogenous assay, add-and-read assay, amenable to HTS format.

Kit Components

For CA-L165

- **Solution I and II:** 5 mL each
- **Component A:** 35 µL
- **Component B:** 120 µL

For CA-L165-10

- **Solution I and II:** 50 mL each
- **Component A:** 350 µL
- **Component B:** 1.2 mL

Storage

Store all in freezer (-20 °C)

Assay Protocol (for one 96-well or 384-well plate)

A) Preparation of Luciferase Assay WORKING SOLUTION

1. Thaw Solution I and Solution II at room temperature (RT, ~22 °C).
2. Thaw components A & B at RT and then keep them on ice. **Note: Do not thaw Component A & B at 37°C.**
3. Prepare Assay Buffer by mixing Solution I (5 mL) and Solution II (5 mL). Transfer Component A (35 µl) and Component B (120 µl) into the 10 ml of Assay Buffer to make luciferase assay WORKING SOLUTION. The amount of WORKING SOLUTION prepared can be modified to suit the experimental design. *Note: Unused Components A and B can be refrozen (at -20 °C). Unused Assay Buffer can be stored at 4 °C to be used for later experiments the same day.*
4. Mix thoroughly by inverting and vortexing for 30 sec. The reconstituted WORKING SOLUTION at RT should be used within 1 hour, or it can be kept at 4°C for no more than 4 hours. In general, only the freshly prepared WORKING SOLUTION is recommended for the experiments.

B) Measurement of Luciferase Activity in Growing Cells

1. Grow cells on opaque white 96 well plates* appropriate for luciferase signal measurement.
*For example: Falcon® 96-well White/Clear Flat Bottom TC-treated Microplate, w/ Lid, Sterile, Item # 353377
2. To generate assay signal, add an equal volume of luciferase assay WORKING SOLUTION to culture media overlaying cells.
3. Incubate the plate at room temperature (~22 °C) for 5 minutes.
4. Read on a luminescence plate reader and record the data.



Label License

The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product for research purposes only by the buyer, where such research does not involve testing, analysis or screening services for any third party in return for compensation on a per test basis. The buyer cannot sell or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise use this product or its components or materials made using this product or its components for Commercial Purposes. Commercial Purposes means any activity by a party for consideration and may include, but is not limited to: (1) use of the product or its components in manufacturing; (2) use of the product or its components to provide a service, information, or data to third parties; (3) use of the product or its components for therapeutic, diagnostic or prophylactic purposes; or (4) resale of the product or its components, whether or not such product or its components are resold for use in research.

